

Scope:

This bulletin covers the new DMSO Filter Support, a product enhancement for the AXP Processing Bagset. (8-5101).

Background:

A new accessory for the AXP device, the DMSO Filter Support, has been introduced as a product enhancement. The use of the accessory is optional and used at the discretion of the end user.

ThermoGenesis Corp has found through Validation and testing that the effectiveness of the DMSO Filter Support will enhance the use of the AXP and minimize filter related issues. The ease of use of the DMSO Filter Support will help stabilize the DMSO line as well as the DMSO filter during centrifugation and aid in the processing of the AXP.

Two additional steps have been added to the loading and unloading of the AXP processing bag set.

ThermoGenesis is providing you with **reusable** DMSO Filter Supports (included with this technical service bulletin). Note: Additional DMSO Filter Supports may be obtained through ThermoGenesis by ordering part no. 0??????????????.

Procedure:

1. New DMSO Filter Support.



2. Connect the DMSO Filter Support on to the tubing, with the notched side against the DMSO filter.



3. Place the DMSO filter into its appropriate place in the AXP with the bottom of the DMSO Filter Support flat against the lower part of the AXP.



4. Proceed with sample processing.
5. Remove the DMSO Filter Support for reuse.

Contact Information:

If you have any questions, contact THERMOGENESIS CORP. Technical Service at 800-783-8357 (U.S. and Canada) or 916-858-5100 (non-U.S./Canada), fax to 916-858-5199, or email at support@thermogenesis.com for assistance.