Scope:

This bulletin covers the introduction of a new accessory for the AXP device, the Component Retaining Clip.

Background:

A new accessory for the AXP device, the Component Retaining Clip, has been introduced as a product enhancement. This clip will help stabilize bag set components stored behind the RBC bag during centrifugation. Two additional steps have been added to the loading and unloading of the processing bag set from the device.


Procedure:

See the attached Instructions for Use (IFU), 37-00-185. After loading the processing bag set into the AXP device, install the Component Retaining Clip over the tubing between Tubing Clamp 2 and Sampling Site C by slipping the clip's hook end over the top edge of the device and snapping the horizontal bar down (see photo in IFU). After centrifugation, the clip can be removed by lifting the horizontal bar up and out of the device (see photo in IFU).

Contact Information:


If you have any questions, contact THERMOGENESIS CORP. Technical Service at 800-783-8357 (U.S. and Canada) or 916-858-5100 (non-U.S./Canada), fax to 916-858-5199, or email at support@thermogenesis.com for assistance.

Component Retaining Clip, IFU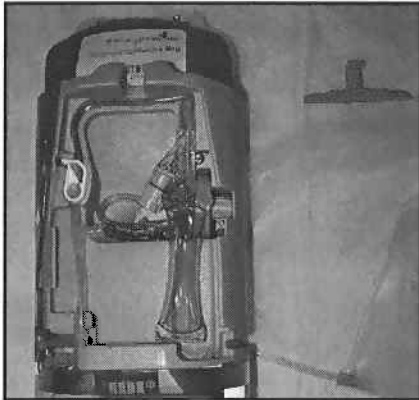

1

ADD CLIP
 Ajouter la fixation
 Aggiungi graffa
 Adicionar clipe
 Lägg till klämma
 Clip hinzufügen
 Clip toevoegen
 Añadir clip

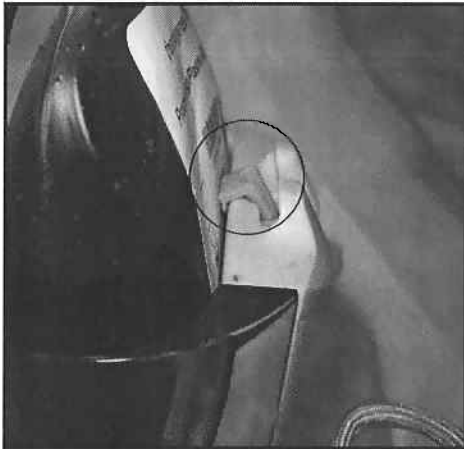2

3

REF 8-5148

INTENDED USE: For use with the AXP Auto Xpress™ Platform

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Fabricado na E.U.A. / In U.S.A. hergestellt
Geproduceerd in U.S.A. / Tillverkad i U.S.A.
ThermoGenesis Corp.
U.S.A